




Tech Tips

tech@rbdinstruments.com

www.rbdinstruments.com

11-155 Power Supply

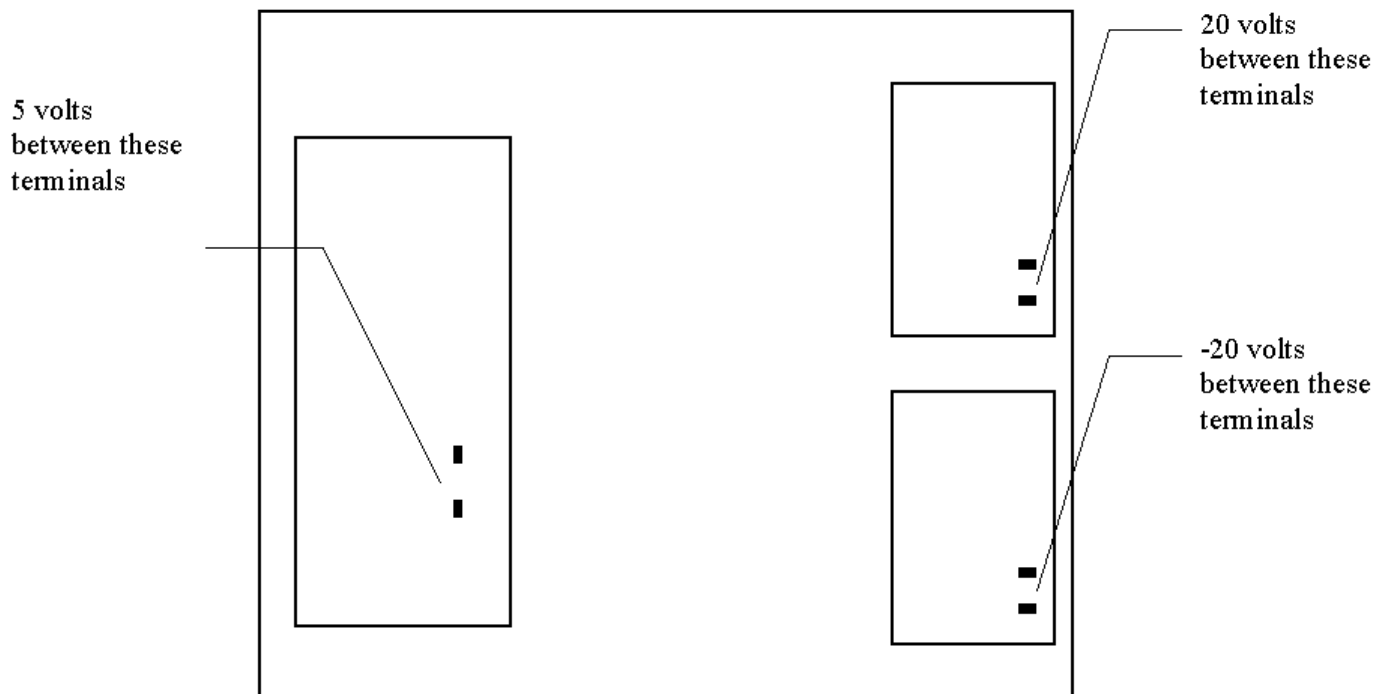

Procedure:

Slide the 11-155 out so that the front of the unit is resting on the floor, and the back of the unit is still in the rack. Remove the top cover.

Turn on the RBD110 main power and measure the voltages as shown in the above drawing. The 5 volt supply usually is between 5 and 5.4 volts DC. The plus and minus 20 volts supplies are usually between 19.5 and 20.5 volts DC. All supplies should have less than 10 mV AC ripple.