



## Tech Tips

[tech@rbdinstruments.com](mailto:tech@rbdinstruments.com)

[www.rbdenter.com](http://www.rbdenter.com)

### **20-150/20-250 Modifications and the RBD147 Interface Unit**

If a system is upgraded to a RBD147 from a PHI PC137A, then the Digital Circuitry Modifications to the 20-150/20-250 need to be removed. The Digital Circuitry Modifications were used on the earlier upgrades, but they cause instability when used in conjunction with the RBD147 interface unit.

All 20-150/250 units no longer have the Digital Circuitry Modification performed, only the video circuitry modification.